

Micro-Fibers Evaluation Utilizing FTIR Microscopy

No. FTIR-2004

■ Introduction

Forensic investigations commonly involve the collection of textile fibers when examining crime scenes. It is vital to characterize and identify those fibers as they can provide extensive information relating to a crime, tying a suspect to a location. Microscopic fibers can present problems in terms of both size and quantity of sample. Fourier transform Infrared spectroscopy (FTIR) presents a nondestructive method to identify those fibers. FTIR can be coupled with microscopy presenting the ability to provide a forensic examiner with both detailed visual microscopic information and simultaneous material identification.

Douglas Deedrick¹ from the FBI stated, "The likelihood of two or more manufacturers duplicating all aspects of the fabric type and color exactly is extremely remote," further exemplifying the importance of fibers at a crime scene and the need to characterize them. Infrared spectroscopy is invaluable for determining the material properties of single fibers.^{2,3} Whether the fibers of interest are synthetic, natural or a hybrid of the two in composition, they will exhibit unique spectral features when measured under infrared light. Fundamental vibrations caused by the absorption of infrared light correspond to structural composition that make up the fibers and dyes commonly used. The advantage of the Shimadzu's compact IRSpirit FTIR spectrophotometer coupled with SurveyIR FTIR microscope to assist the forensic examiner in characterizing microscopic fibers is presented in this application note.

■ Results and Discussion

The IRSpirit FTIR spectrophotometer and SurveyIR™ FTIR microscope from Redwave Technologies provide a simplified field-deployable system for fiber analysis and identification. Since the microscope utilizes the standard detector of the FTIR, the need for liquid nitrogen for cooling the MCT detector may be eliminated.

The live image feed from the software allows for simultaneous viewing and spectral data collection. The SurveyIR microscope can switch between reflectance and transmittance collection modes as well as a clip-on view-through attenuated total reflection (ATR) accessory.

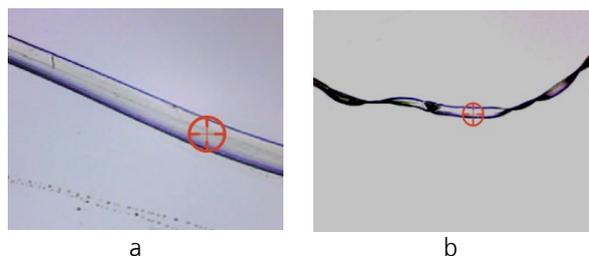


Figure 1: SurveyIR images captured of two synthetic fibers, polyester (a) and polyamide (b). Red crosshairs indicate the sampling position at an aperture size of 60µm.

The images in figures 1 and 3 represent fibers extracted from woven fabric. The fibers were cut into small lengths (1-3mm), rolled flat via roller knife on plain glass microscope slides, and then transferred to Low-E glass microscope slides for analysis. The left image of polyamide (Fig 1, a) shows a transparent fiber about 60µm in diameter. The right image of polyester (Fig 1, b) is slightly less than 60µm wide. Due to the sample preparation technique used, structural appearance of these two fibers is very similar as both fibers appear to be similar in size and transparency; however, their corresponding IR spectra clearly represent two different materials.

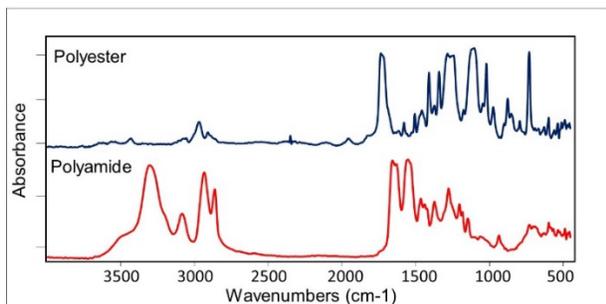


Figure 2: Overlaid spectra of polyester (blue) and polyamide (red) fibers.

FTIR spectra were collected by reflectance at 8cm^{-1} resolution and coadded 64 scans. The bottom spectrum of Figure 2, of a polyamide, exhibits spectral features consistent with nylon 6,6, poly(hexamethylene apidamide). The top spectrum of Figure 2 is consistent with polyethylene terephthalate (PET). Both fibers can be differentiated utilizing FTIR spectroscopy and identified using spectral library searching methods.

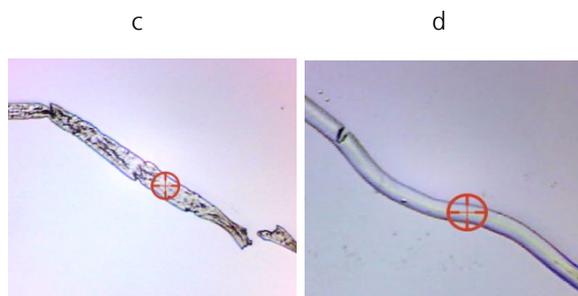


Figure 4: SurveyIR images captured of two natural fibers, cotton (left, c) and silk (right, d). Red crosshairs indicate sampling are with both apertures set to $60\mu\text{m}$.

Cotton and silk are well known textiles and some of the oldest known. Cotton has been found in archeological sites dating back as early as 5000-6000 BC. Likewise, silk has history dating back to ancient Chinese civilizations, usually in the hands of the emperors or elite members of society. Both materials have very different textures, as cotton is known to be soft and fluffy where silk is very smooth and lustrous. Unlike the synthetic fibers, the images in Fig. 3 exhibit apparent differences as silk (d) appears very smooth compared to cotton (c) which has small inconsistencies. The composition of the two is completely natural; however, cotton is comprised of almost completely of cellulose and silk, protein.

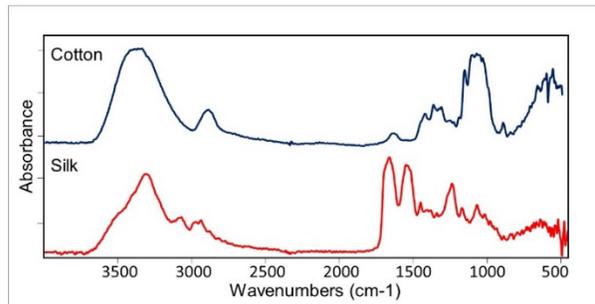


Figure 3: Overlaid spectra of cotton (blue) and silk (red) fibers.

The IR spectra seen in Fig. 3 of silk (red) has distinct peaks associated with N-H structural units around 3300cm^{-1} and amide bands I and II at 1661 and 1532cm^{-1} , respectively. Cotton (Fig. 3, blue) on the other hand has an absence of amide bands I and II (peaks 1661 and 1532cm^{-1}) and a broadened peak at 3380cm^{-1} due to extensive hydrogen bonding in the cellulose structure. The broaden peak at 1085cm^{-1} (blue) is associated with cellulose structure as a combination of C-C, C-OH, C-O-C stretches. The ability of FTIR to differentiate the two natural fibers is apparent and can be referenced against a library for identification purposes.

■ Conclusion

FTIR microscopy utilizing IRspirit FTIR spectrophotometer and SurveyIR FTIR microscope (Figure 5) proves to be a useful tool in analyzing and identifying small fibers. Single fiber analysis can be accomplished without sample destruction and reference images captured for later presentation and sample position validation. The resulting spectra from both natural and synthetic fibers were identified utilizing library searches and further verified by key characteristic fundamental IR vibrations.

The uniqueness of IR vibrational spectroscopy is proven in both synthetic and natural fibers analyzed providing certainty for identification purposes. To the forensic examiner, certainty is imperative for evidence that is to be used in the court. The ability of SurveyIR to switch between different sampling methods (ATR, Transmittance, Reflectance) presents an effective, easy to use toolset identify and differentiate materials with high confidence during forensic analysis.



Figure 5: IR Spirit FTIR Spectrometer with SurveyIR Microscope Accessory with Diamond ATR attachment.

■ Acknowledgement

This document was created as a result of collaborative work with Redwave Technologies (formerly Czteck Inc.). Permission to modify Redwave Technologies' original document: "Evaluation of Micro-Fibers Utilizing Microspectroscopy", was obtained. We are grateful to them for accommodating our request.

■ References

1. Deedrick, Douglass W. "Hairs, Fibers, Crime, and Evidence Part 2: Fiber Evidence," *Forensic Sci. Comm.*, **2**, 3 (2000).
2. ASTM E2224 - 10 , "Standard Guide for Forensic Analysis of Fibers by Infrared Spectroscopy," ASTM International, 100 Barr Harbor Drive, PO Box C700, West Conshohocken, PA.
3. "A FORENSIC FIBER EXAMINER TRAINING PROGRAM," Scientific Working Group for Materials Analysis (SWGMA), <https://archives.fbi.gov/>



SHIMADZU Corporation
www.shimadzu.com/an/

SHIMADZU SCIENTIFIC INSTRUMENTS
7102 Riverwood Drive, Columbia, MD 21046, USA
Phone: 800-477-1227/410-381-1227, Fax: 410-381-1222
URL: www.ssi.shimadzu.com

First Edition: August 2020

For Research Use Only. Not for use in diagnostic procedure.
This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. Shimadzu disclaims any proprietary interest in trademarks and trade names used in this publication other than its own. See <http://www.shimadzu.com/about/trademarks/index.html> for details.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and