

Application News

Material Testing System

MCT

Compression Test for Structural Materials of Lithium-Ion Batteries by MCT

No. SCA_300_002

Introduction

Since lithium-ion batteries are light and small, they are used in a wide variety of products, from mobile electronic devices such as cellular phones and notebook PCs to electric cars and hybrid cars. Their inner structural materials are subjected to external force during production processes and to pressure during use. Therefore, evaluation of strength of each structural material is important to maintain consistent quality. A strength measurement was performed on thin or minute materials among various structural materials of lithiumion batteries. Separators are usually evaluated by a tensile test or penetration test. A compression test is also important to evaluate them because they are compressed in some processes. Active materials of approximately $10 \,\mu$ m in size located near the electrode need to have a certain compression strength so they will not be destroyed during the coating process. Below are the results of compression tests performed on these materials using the MCT-211 Series Micro Compression Testing Machine.



Fig. 1 External View of the MCT-211 Series (with the Side Observation Kit Mounted)



Fig. 2 Structure of Lithium-Ion Battery

■Compression Test on Separator of Lithium-Ion Battery

Table 1 shows the three types of specimens used for the measurement. Table 2 shows accessories used in the test and test conditions. Fig. 3 shows the conceptual diagram of the measurement. Table 3 shows the results of the compression tests on the three types of specimens. The specimens were evaluated by a compression rate where the same test force was applied. The results clearly show the difference among the three types. Fig. 4 is a graph indicating the test force - displacement relationship of each specimen. The inflexion point of specimen 2 is at around 10 mN (pressure of approximately 5 MPa), indicating that applying too much compression pressure causes plastic deformation to the separator.

Table 1	Specimens
---------	-----------

1) Specimen Name	Separator		
2) Specimen Number	1	2	3
3) Thickness	20 µm	20 µm	10 µm

	Conditione
1) Upper Indenter	Flat indenter (with a diamond tip), 50 μm dia.
2) Test Mode	Load-unload test
3) Test Force (mN)	50
4) Loading Rate (mN/sec)	2.2
5) Holding Time (sec)	0
6) Test Method	A thin layer of liquid glue was applied to a glass plate, the separator was bonded to it, and a compression test was performed using the upper indenter. (See Fig. 2.)

Table 2 Test Conditions



Fig. 3 Conceptual Diagram of Measurement

Table 3 Test Results

Specimen Name	Specimen Number	Maximum Force (mN)	Compression Variation (µm)	Compression Rate (%)
	1	49.9	3.651	18.3
Separator	2	49.9	3.371	16.9
	3	50.0	1.038	10.4

Note: The compression rate was calculated by the following calculation formula. Compression rate $(\%) = (\text{compression volume})/(\text{thickness}) \times 100 (\%)$



Fig. 4 Results of Compression Test

■Compression Test on Active Materials

A compression test was performed on two types of active material particles for positive electrode of lithium-ion battery. Table 4 shows test conditions and Fig. 5 shows an image of the test (compressed part). Measurement was performed ten times for each specimen. Then, average values were selected as a representative value for each specimen. (See Table 5 and Fig. 6.) The results clearly indicate the difference in the strength between the two active materials and lithium cobalt oxide (LiCoO2) was confirmed to have the higher strength. As shown above, the MCT-211 Series Micro Compression Testing Machine enables accurate and efficient evaluation of compression characteristics of thin or minute materials used inside lithium-ion battery.

1) Upper Indenter	Flat indenter (with a	
	diamond tip),	
	50 µm dia.	
2) Test Mode	Compression test	
3) Test Force (mN)	50	
4) Loading Rate	2.2	
(mN/sec)		
5) Holding Time (sec)	0	
6) Test Method	A very small amount of	
	each specimen was spread	
	on the lower compression	
	plate and a compression	
	test was performed on each	
	single particle. (See Fig. 5.)	

Table 4 Test Conditions



Fig. 5 Conceptual Diagram of the Compressed Part

Table 5 Test Results

Specimen Name	Fracture Force (mN)	Particle Diameter (µm)	Strength (MPa)
LiMn ₂ O ₄	1.67	13.0	7.79
LiCoO ₂	16.23	13.3	72.75

Note: Fracture strength was calculated by the following calculation formula

(Based on JIS R1639-5*1).

 $Cs = 2.48 P/\pi d2$

Cs: Strength (MPa), P: Fracture force (N), d: Particle diameter (mm)

*1: Test methods of properties of fi ne ceramic granules Part 5: Compressive strength of a single granule



Fig. 6 Force - Displacement Curves



Shimadzu Europa GmbH www.shimadzu.eu

For Research Use Only. Not for use in diagnostic procedures. The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to be the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.