


LC-10ADvp

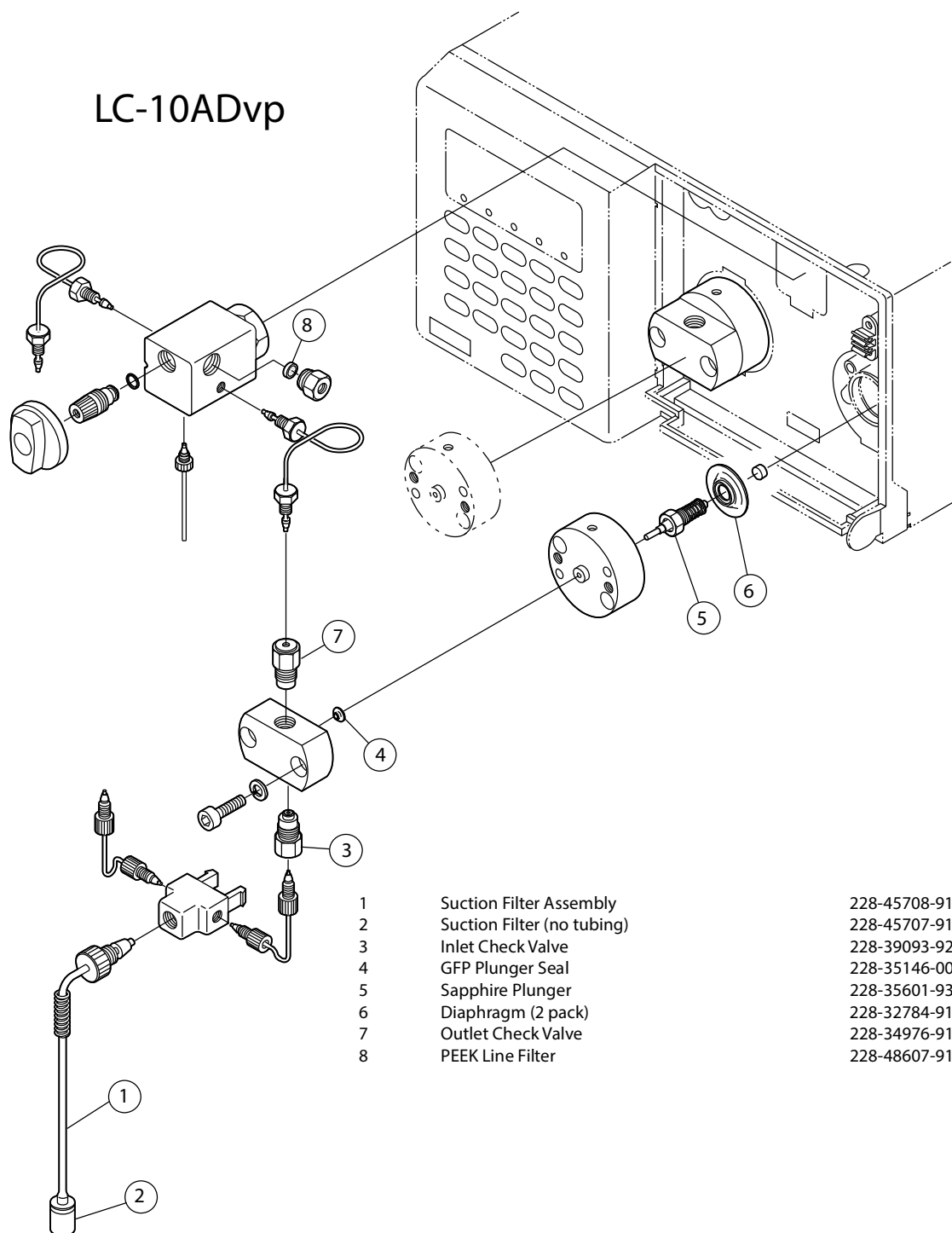

- | | | |
|---|----------------------------|--------------|
| 1 | Suction Filter Assembly | 228-45708-91 |
| 2 | Suction Filter (no tubing) | 228-45707-91 |
| 3 | Inlet Check Valve | 228-39093-92 |
| 4 | GFP Plunger Seal | 228-35146-00 |
| 5 | Sapphire Plunger | 228-35601-93 |
| 6 | Diaphragm (2 pack) | 228-32784-91 |
| 7 | Outlet Check Valve | 228-34976-91 |
| 8 | PEEK Line Filter | 228-48607-91 |